LCV TEST KİTİ

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LCV TEST KIT

Leuco- crystal violet (LCV) is a chemistry widely used in laboratory and field studies to show the presence of blood. The leuco-crystal violet (LCW) reacts with the heme group to give a violet color.

With this test kit, blood stains can be determined easily and quickly using the LCV method.

CONTENT OF TEST KIT	
Alcohol Reagent	30 mL
LCV Reagent	30 mL
Hydrogen Peroxide Reagent	30 mL
Distilled Water –15 mL	
Swab – 50 Pcs	
Filter Paper – 50 Pcs	
Example of known Blood Stain - 1 Pc	

LCV TEST METHOD

❏ Sampling:

A drop of **Alcohol Reagent** is placed on a cotton swab or filter paper and then rubbed onto a stained surface.

❏ PROCEDURE:

Add 1 drop of **LCV Reagent**, and 1 drop of **Hydrogen Peroxide Reagent** onto the sampled cotton swab or filter paper.

❏ RESULT:

- **Positive Test:** If the color of the cotton swab or filter paper turns blue - purple immediately, the result is interpreted as **POSITIVE TEST**. So the sample is blood.
- **Negative Test:** If the color of the cotton swab or filter paper lasts for 1 minute or longer, the result is interpreted as **NEGATIVE TEST**. So the sample is not blood.
- You can observe a positive result by experimenting with a known blood stain sample before performing the experiment.